



Glass Plating Beads (Sterilized)

FYO001-100G

Glass Plating Beads (Sterilized) provide a easy way to perform a routine procedure plating bacteria or yeast. The beads can be recovered by washing with 70% ethanol and sterilized by autoclave and dried in oven. Each package contains enough beads for 100 to 200 plates.

Protocol for plating bacteria and yeast

Pipet the cells on the plate, add 5 to 10 beads, and move the plate back and forth a few time. The beads' rolling action gently spreads the cells evenly and completely over the agar surface. Entire stacks of plating, pour the beads off the plate and it is ready for the incubation.